Autoclavable Micro Pipette

Nichipet® Premium L




Innovation in "Fatigue-Free" & "Comfort"



What's new!





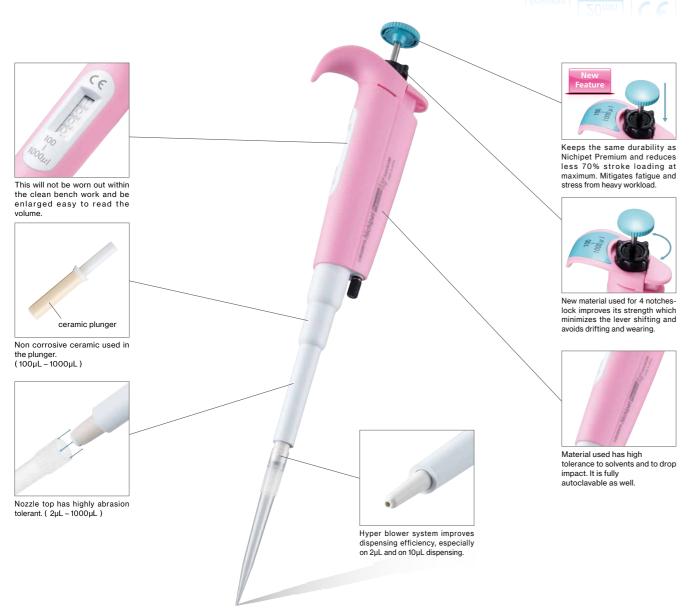

- ■The durability still is same as Nichipet Premium has. Stroke loading of Nichipet Premium LT becomes less 70% at maximum than before.
- ■Hand temperature doesn't permeate through the body of the pipette which works to avoid affecting volume measurement. (Patented)
- *All features of Nichipet Premium are available on Nichipet Premium LT.
- Since the volume indicator is made of glass, ultraviolet germicidal of the clean bench will not make it cloud.
- From prolonged irradiation, the surface layer of plastic may deteriorate or discolor, but its performance is not affected.

Cat. No	Volume Setting	Volume range(µL)	Increments(µL)	Usable tips			
00-NLT-2	Digital Setting	0.1 – 2	0.002	BMT2-UT/UTWR/UTWRB、FUT/FUTRB			
00-NLT-10		0.5 – 10	0.01	BMT2-SSW/SSWR/SSWRB、FSS/FSSRB			
00-NLT-20		2-20	0.02	BMT2-SG/SGR/SGRB、FSG/FSGRB			
00-NLT-100		10 – 100	0.1	BMT2-SG/SGR/SGRB、FSE/FSERB			
00-NLT-200		20 – 200	0.2	BMT2-SG/SGR/SGRB、FAG/FAGRB			
00-NLT-1000		100 – 1000	1	BMT2-LG/LGR/LGRB、FLG/FLGRB			

^{*}See p.21-22 for tip selection.

Accuracy and Precision

Cat. No	Volume(µL)	Accuracy(%)	Precision(%)	Cat. No	Volume(µL)	Accuracy(%)	Precision(%)
00-NLT-2	0.2	±12.0*	≦6.0*	00-NLT-100	10	±2.0	≦1.0
	1	±5.0	≦2.5		50	±1.0	≦0.3
	2	±3.0	≦1.0		100	±0.8	≦0.3
00-NLT-10	1	±4.0*	≦3.0 *	00-NLT-200	20	±1.0	≦0.5
	5	±1.0	≦1.0		100	±0.8	≦0.3
	10	±1.0	≦0.5		200	±0.8	≦0.2
00-NLT-20	2	±5.0	≦3.0	00-NLT-1000	100	±1.0	≦0.5
	10	±1.0	≦1.0		500	±0.8	≦0.3
	20	±1.0	≦0.4		1000	±0.7	≦0.2

^{*}Each accuracy and precision are the values which come from measuring with our compatible tip.
*The AC and CV of NLT-2 below 0.2 \(\mu \) L and NLT-10 below 1 \(\mu \) L depend on the sampling skill and the sampling condition where pipettes are used.

Nichiryo's quality starter packages

Nichipet Premium LT S-PAC

- Package type of 3 selected Pipettes with racked -Tips.
- Ready-to-start packages with 3 sizes of pipettes to perform high AC and CV pipetting.
- ■3 packages are available to cover wide range of volume.





Racked Tip 3 or 4 racks







Nichiryo Tube Opener (code:00-NTO)

"Easier" and "safer" to open tube lid with holding pipette



- ■Attached on grip end, 0.5, 1.5 mL tube and screw cap can be opened easily.
- *Compatible with Nichipet Premium LT ($2\text{-}1000\mu L$) and Nichipet EX ${\rm I\!I}\,$ ($2\text{-}1000\mu L$)

Nichipet Premium LT S-PAC

Description	Cat. No	Content
Nichipet Premium LT S-PAC A	00-SP-AL	NLT-10、NLT-1000, NLT-1000 BMT2-SSWR(0.5~10 μ L) 、BMT2-SGR(2~200 μ L) 、BMT2-LGR(100~1000 μ L) 、NTO
Nichipet Premium LT S-PAC B	00-SP-BL	NLT-20 NLT-200 NLT-1000 BMT2-SGR(2 \sim 200 μ L) ×2 BMT2-LGR(100 \sim 1000 μ L) NTO
Nichipet Premium LT S-PAC H	00-SP-HL	NLT-2、NLT-10、NLT-100 BMT2-UTWR(0.1~10 μ L) \times 2、BMT2-SSWR(0.5~10 μ L) 、BMT2-SGR(2~200 μ L) 、NTO