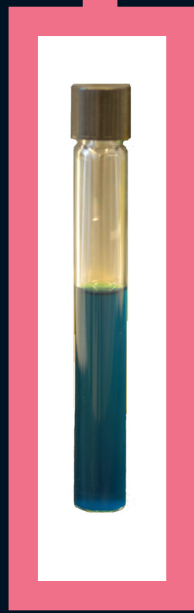
## RambaQUICK™ Salmonella Method

Broth enrichment  
RambaQUICK™ Salmonella

+

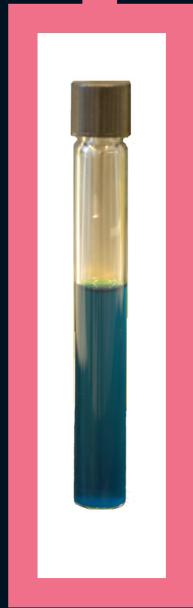
CHROMagar™ Salmonella Plus  
Culture media


Click below:

# ● RambaQUICK™ Salmonella Method

+

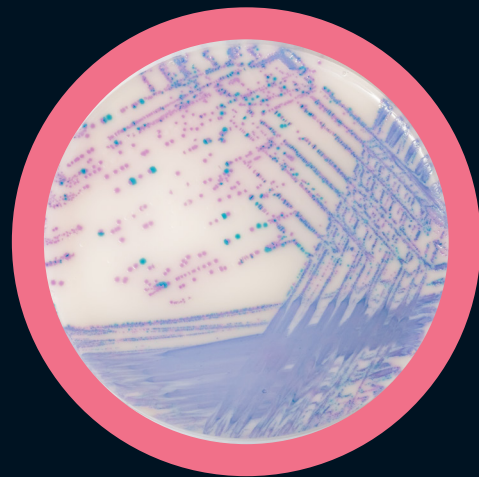Broth enrichment  
RambaQUICK™ Salmonella

CHROMagar™ Salmonella Plus  
Culture media


For Rapid detection of *Salmonella* species  
including *S. typhi*, *S. paratyphi* and  
lactose positive *Salmonella*

# RambaQUICK™ Salmonella Method

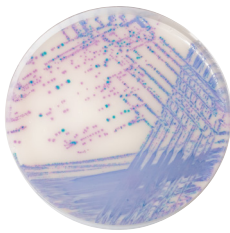www.CHROMagar.com

Broth enrichment  
RambaQUICK™ Salmonella

CHROMagar™ Salmonella Plus  
Culture media

For Rapid detection of *Salmonella* species including *S. typhi*, *S. paratyphi* and lactose positive strains in food samples

## Background

From the CDC report Trends in Foodborne Illness, 1996-2010:

« This year's report (released in CDC's Morbidity and Mortality Weekly Report [MMWR]) summarizes 2010 data from FoodNet and provides the best measure of trends in foodborne disease from 1996–2010. It confirms that *Salmonella* infection has not declined in 15 years,... »  
One of the report recommendations to reduce the number of *Salmonella* infections is "Enhancing laboratory testing and disease reporting to more quickly identify outbreaks and their causes »

## RambaQUICK™ Salmonella Method: Rapid and Simple !

The RambaQUICK™ Salmonella method was designed with the aim of **simplifying the test procedure while maintaining a high level of sensitivity within a conventional culture methodology.**

### 1 TWO IN ONE ENRICHMENT STEP

The RambaQUICK™ selective broth is a combination of the best functions found in each of both conventional enrichments, the RVS and the MKTTn broths. After the revivification phase in BPW, the *Salmonella* replicates at the exponential growth phase in the optimised RambaQUICK™ Salmonella broth, which offers not only a highly nutrient environment but also inhibits the growth of competitive flora.

### 2 FAST METHOD

- negative in 2 days.
- **positive** in 3 days.

### 3 SIMPLE

- one selective enrichment.
- one culture plate.

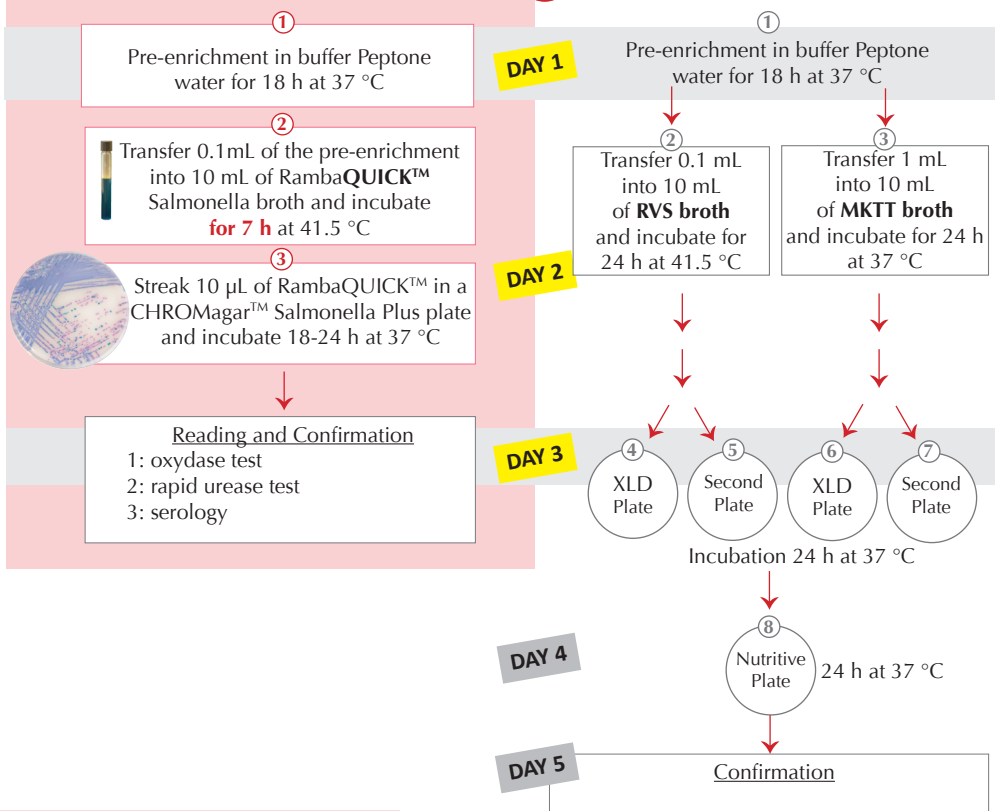
### 4 MEETS ISO 6579 REQUIREMENTS

by detecting lactose positive *Salmonella* in intense colour on CHROMagar™ Salmonella Plus.

- *Salmonella* including *S. typhi*, *S. paratyphi* A and lactose positive *Salmonella* → Mauve

## RambaQUICK™ Salmonella Method

VS ISO standard method ISO 6579

UP TO 11 DAYS  
FOR POSITIVE SAMPLES

## Method Description

RambaQUICK™ Enrichment Broth Powder	Total .....	34.2 g/L	
	Salts .....	8.6	
+	Peptone .....	4.5	
	Selective mix .....	21.1	
	Powder Storage at 15/30°C - pH: 5.6 +/-0.2 Shelf Life: 2 years		
	CHROMagar™ Salmonella Plus Culture Media Powder	Powder Base : Total .....	32.8 g/L
	Agar .....	15.0	
	Peptone and Yeast extract ...	8.0	
	Salts .....	8.5	
	Chromogenic mix .....	1.3	
	Storage at 15/30 °C - pH: 7.5 ± 0.2 Shelf Life: 3 years		
	Supplements	1* : Liquid form .....	6.0
		Storage at 15/30 °C Shelf Life: 5 years	
		In order to obtain a white opaque background:	
		2 <sup>nd</sup> : White Opaque Supplement (Powder):	1.0
		Storage at 15/30 °C Shelf Life: 3 years	

Scientific Publications on this product: available on [www.CHROMagar.com](http://www.CHROMagar.com)

Please read carefully the instructions for use (IFU document) available on [www.CHROMagar.com](http://www.CHROMagar.com)

## Order References

Please use these product references when contacting your local distributor:

RambaQUICK™ Salmonella Enrichment broth powder in 1 000 mL pack (100 tests of 10 mL): SQ001  
CHROMagar™ Salmonella Plus culture media powder in 5 000 mL pack (250 plates of 20 mL): SA162

RambaQUICK™ Salmonella Method

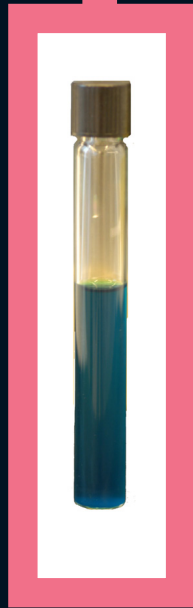
to prepare 5 000 mL of media (250 tests): 3 packs of SQ001 + 1 pack of SA162

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